


FID/NPD Adaptor Fitting for Agilent GCs

(cat.# 20884)

Installation Instructions

1. Cool the GC and remove the old detector fitting.
2. Assemble the fitting as shown and screw tightly into the GC.
3. Insert the column 28mm from the back of the nut. Alternatively, install the column until it touches the jet restriction, then pull back approximately 1mm to allow for forward movement when the ferrule is compressed.
4. Leak check with the Restek Leak Detector (cat.# 22839).

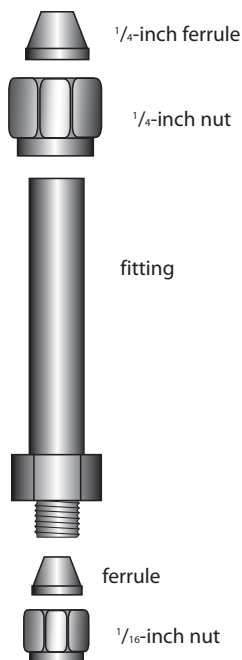

Restek trademarks: Restek logo

***Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4
(or your Restek representative) if you have any questions about
this product or any other Restek product.***



#203-03 [005]
Rev: 07/2008

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823
phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • www.restek.com
Restek France • phone: 33 (0)1 60 78 32 10 • e-mail: restek@restekfrance.fr
Restek Ireland • phone: 44 2890 814576 • e-mail: restekeurope@aol.com
Thames Restek U.K. LTD • phone: 44 1494 563377 • e-mail: sales@thamesrestek.co.uk
Restek GmbH • phone: +49 (0) 6172 2797 0 • e-mail: info@restekgmbh.de