


Restek MSD Conversion Fitting

(cat.# 21314)

This MSD Conversion Fitting is used to convert your source fittings for use with standard nuts and ferrules. This design uses Vespel®/graphite ferrules and inexpensive 1/16-inch nuts instead of the costly instrument manufacturers capillary nuts. The kit includes the MSD conversion fittings and aluminum sealing rings.

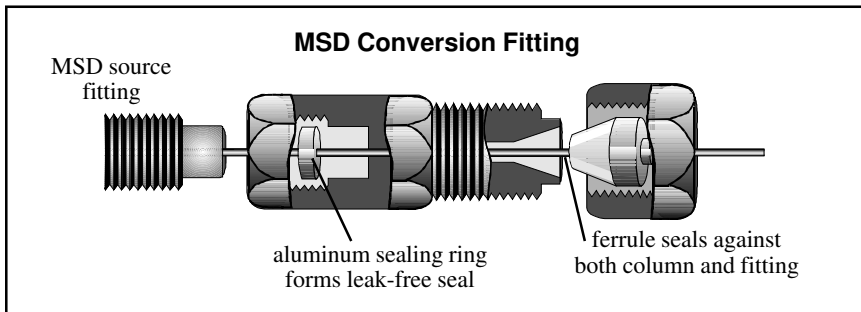

Installation Instructions

1. Install aluminum sealing ring into the MSD adaptor fitting.
2. Screw the MSD Conversion Fitting onto the MSD source. Be sure to use a wrench on the MSD source while tightening down the fitting to avoid damage to the MSD source.
3. After installing the MSD Conversion Fitting, place a no-hole ferrule in the 1/16-inch nut and check for leaks by monitoring mass 32 or 28.
4. After the system has been certified leak-free, use the appropriate Vespel® ferrule to seal the column. Always place a wrench on the MSD Conversion Fitting when tightening or loosening.

Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your local Restek representative) if you have any questions about this product or any other Restek product.



#203-03 [013]
Rev. date: 3/01

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: dmannusrestek@compuserve.com