


Finger-Tight Column Nuts

For use with Agilent®-style ferrules - cat.# 21040

Recommended Installation Distance:

Restek recommends an installation distance of 26–28 mm from the back of the nut to the end of the fused silica column.

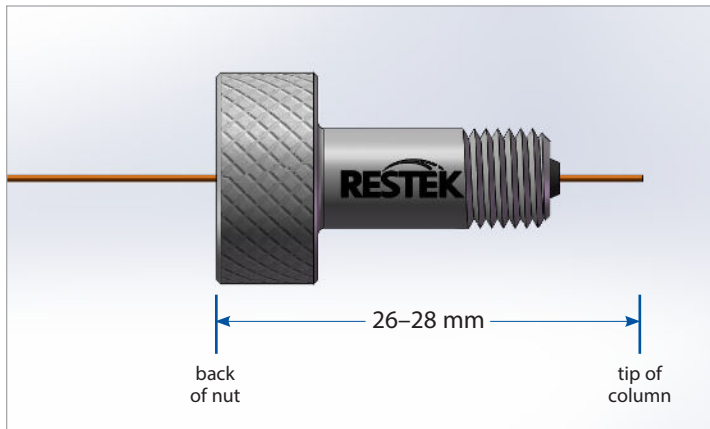

Questions about this or any other Restek® product? Contact us or your local Restek® representative (www.restek.com/contact-us).

Restek® patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.) Other trademarks in Restek® literature or on its website are the property of their respective owners. Restek® registered trademarks are registered in the U.S. and may also be registered in other countries.

© 2014 Restek Corporation. All rights reserved.
Printed in the U.S.A.

#203-03-024 Rev. date: 10/14



RESTEK®
www.restek.com