


ECD Vent Trap Kit

(Includes ECD Vent Trap and 9-inch piece of copper tubing)

cat.# 22017

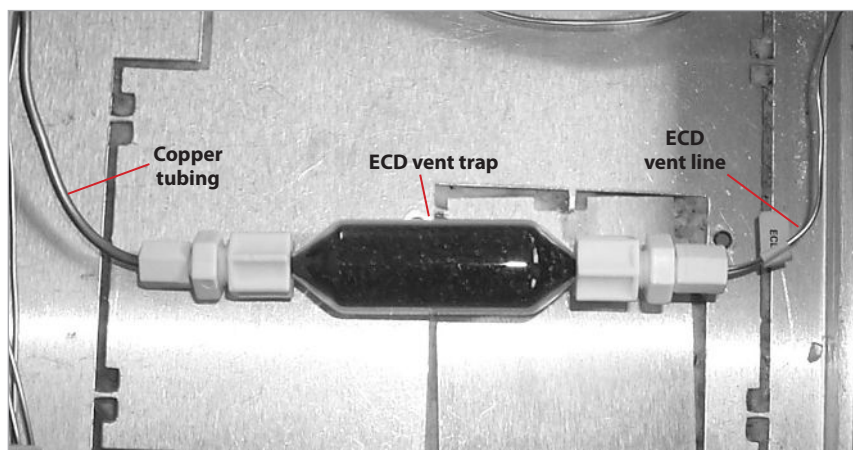

1. Connect the 9-inch piece of copper tubing to one end of the ECD vent trap.
2. Connect the other end of the ECD vent trap to the existing ECD vent line.
3. Position the ECD vent trap in the GC as desired. Avoid placing the trap in locations where it could be broken.

Questions about this or any other Restek product?
Contact us or your local Restek representative (www.restek.com/contact-us).

Restek patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.) Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.

© 2024 Restek Corporation. All rights reserved. Printed in the U.S.A.

www.restek.com

#204-08-009 Rev. date: 12/24



RESTEK
Pure Chromatography