

On-Column Liner for Agilent & CIS4 PTV Inlets

cat.# 23430, 24977

Use **only** 0.53 mm columns for Press-Tight connection in liner


- A 0.53 mm analytical column directly installed in the liner, or a 0.53 mm retention gap connected to an analytical column (0.53 mm or less) is necessary.
- Smaller ID columns will not allow the syringe needle to enter the head of the column.

Use **only** tapered 23-26s gauge needles with a cone-type needle style

- Larger-sized needles will not enter the 0.53 mm column.
- Cone-type needle style will deliver the sample directly into the column.

With autosamplers, use **only** slow injection speed option

- This will minimize the likelihood of syringe damage.

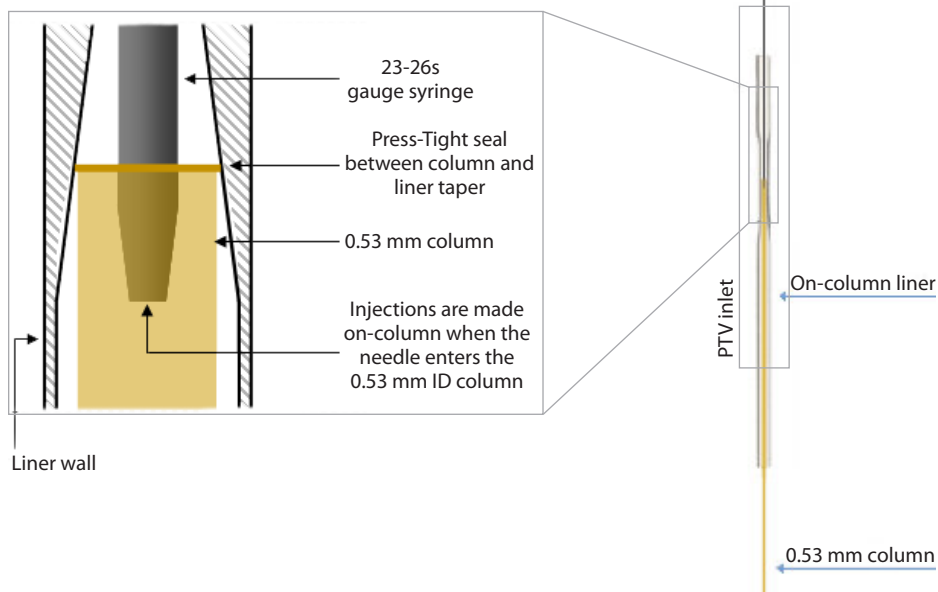

Questions about this or any other Restek product? Contact us or your local Restek representative (www.restek.com/contact-us).

Restek patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.) Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.

© 2020 Restek Corporation. All rights reserved. Printed in the U.S.A.

www.restek.com

#205-02-001 Rev. date: 09/20



RESTEK
Pure Chromatography