

SS Secure-Fit Fittings

Single ended: cat.#s 25181–25193

Double ended: cat.#s 25208—25216

Secure-Fit connectors are designed to ensure a leak-free seal and eliminate dead volume with any standard $1/16"$ - $10/32$ internal HPLC fitting connection port.

Instructions for Use:

- Push the Secure-Fit fitting into the port until the threads engage.
- Turn the Secure-Fit fitting until it is finger-tight. Secure-Fit fittings seal easily so only a light tightening is required to obtain a leak-free seal. The tubing penetration depth is automatically set.
- For double ended fittings, repeat the above steps.
- For single ended fittings, be sure to make a good zero dead volume connection at the other end of the tube with standard nuts and ferrules.



***Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4
(or your local Restek representative) if you have any questions
about this product or any other Restek product.***



ISO 9001 #730-20 [001]
Rev. date: 12/28/01

Restek USA: 110 Benner Circle, Bellefonte, PA 16823
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; www.restekcorp.com
Restek GmbH: phone: (49) 06172 2797 0 • email: RESTEK-GMBH@t-online.de
Restek France: phone: (33) 01 60 78 32 10 • email: restekfr@club-internet.fr
Thames Restek UK Ltd.: phone: (44) 01753 624111 • email: Sales@Thamesrestek.co.uk
Thames Restek Scotland Ltd.: phone: 0870 241 1247 • email: Sales@Thamesrestek.co.uk
Restek Ireland: phone: (44)2 890 814 576 • email: restekeurope@cs.com