

# PEEK® Secure-Fit Fittings

**Single ended:** cat.#s 25217–25220, 25230–25237


**Double ended:** cat.#s 25221—25229

---

Secure-Fit connectors are designed to ensure a leak-free seal and eliminate dead volume with any standard  $1/16"$ - $10/32$  internal HPLC fitting connection port.

## Instructions for Use:

- Push the Secure-Fit fitting into the port until the threads engage.
- Turn the Secure-Fit fitting until it is finger-tight. Secure-Fit fittings seal easily so only a light tightening is required to obtain a leak-free seal. The tubing penetration depth is automatically set.
- For double ended fittings, repeat the above steps.
- For single ended fittings, be sure to make a good zero dead volume connection at the other end of the tube with standard nuts and ferrules.

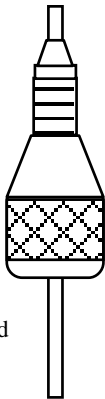

***Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4  
(or your local Restek representative) if you have any questions  
about this product or any other Restek product.***

---



ISO 9001 #730-20 [002]  
Rev. date: 12/28/01

**Restek USA:** 110 Benner Circle, Bellefonte, PA 16823  
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; [www.restekcorp.com](http://www.restekcorp.com)  
**Restek GmbH:** phone: (49) 06172 2797 0 • email: [RESTEK-GMBH@t-online.de](mailto:RESTEK-GMBH@t-online.de)  
**Restek France:** phone: (33) 01 60 78 32 10 • email: [restekfr@club-internet.fr](mailto:restekfr@club-internet.fr)  
**Thames Restek UK Ltd.:** phone: (44) 01753 624111 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)  
**Thames Restek Scotland Ltd.:** phone: 0870 241 1247 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)  
**Restek Ireland:** phone: (44)2 890 814 576 • email: [restekeurope@cs.com](mailto:restekeurope@cs.com)