

# HPLC Fitting Removal Tool

Part #25325

---

## Instructions for Use:

- Drill an opening approximately 1/4" deep in the center of the PEEK fitting by turning the narrow drill bit in a clockwise direction. To prevent damage to the head of the column, do not drill deeper than 1/4".
- Once the opening in the fitting has been created, remove the drill bit.
- Insert the extraction tool into the opening. The extraction tool has a tapered bit and reverse threading.
- Turn the extraction tool in a counterclockwise direction. The normal direction used when **removing** a PEEK fitting. The tool should grab the PEEK fitting and begin to unscrew the fitting from the column.

***Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4  
(or your local Restek representative) if you have any questions  
about this product or any other Restek product.***

---



#730-60 [001]  
Rev. date: 3/6/02

**Restek USA:** 110 Benner Circle, Bellefonte, PA 16823  
phone: (800) 356-1688 or (814) 353-1300 • fax: (814) 353-1309; [www.restekcorp.com](http://www.restekcorp.com)  
**Restek GmbH:** phone: (49) 06172 2797 0 • email: [RESTEK-GMBH@t-online.de](mailto:RESTEK-GMBH@t-online.de)  
**Restek France:** phone: (33) 01 60 78 32 10 • email: [restekfr@club-internet.fr](mailto:restekfr@club-internet.fr)  
**Thames Restek UK Ltd.:** phone: (44) 01753 624111 • email: [Sales@Thamesrestek.co.uk](mailto:Sales@Thamesrestek.co.uk)  
**Restek Ireland:** phone: (44)2 890 814 576 • email: [restekeurope@cs.com](mailto:restekeurope@cs.com)