

Outlet Cap and Gold Seal Assembly Tool

cat.# 24989

Use Outlet Cap cat.# 25139 • Use Gold Disk Seal cat.# 25467



Outlet Cap and Gold Seal Assembly Tool
for Agilent 1050 & 1100 HPLC Systems



Figure 1



Figure 2



Figure 3

Easily install the gold seal into the outlet cap.

1. Put Outlet Cap on male part of Assembly Tool (Figure 1).
2. Push the Assembly Tool together (Figure 2).
3. Hold onto Outlet Cap and pull Assembly Tool apart (Figure 3).
4. Put Gold Seal over pin on male part of Assembly Tool (Figure 4).
5. Push the Assembly Tool together and press the Gold Seal in the Outlet Cap (Figure 5).
6. Pull the Assembly Tool apart and remove assembled Outlet Cap and Gold Seal (Figure 6).



Figure 4



Figure 5



Figure 6

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative)
if you have any questions about this, or any other, Restek product.**



#730-80-005
Rev. date: 02/07

Restek U.S. • 110 Benner Circle • Bellefonte, PA 16823

phone: 814-353-1300 or 800-356-1688 • fax: 814-353-1309 • www.restek.com

Restek France • phone: 33 (0)1 60 78 32 10 • fax: 33 (0)1 60 78 70 90 • e-mail: restek@restekfrance.fr

Restek Ireland • phone: 44 2890 814576 • fax: 44 2890 814576 • e-mail: restekeurope@aol.com

Thames Restek U.K. LTD • phone: 44 1494 563377 • fax: 44 1494 564990 • e-mail: sales@thamesrestek.co.uk

Restek GmbH • phone: +49 (0) 6172 2797 0 • fax: +49 (0) 6172 2797 77 • e-mail: info@restekgmbh.de