

2-in-1 Filter/O-Ring Insertion Tool Kit


for ASE 100/150/200/300/350 Extraction Cells

cat.# 26181–26183

Inserting a filter (ASE 100/150/200/300/350 systems), using Filter Insertion Attachments on 2-in-1 Filter/O-Ring Insertion Tool Handle.



Screw the appropriate size attachment onto the end of the 2-in-1 Filter/O-Ring Insertion Tool Handle.

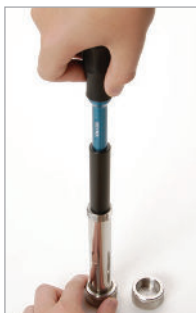

Place a filter at the top of the extraction cell.



Push the filter to the bottom of the extraction cell.



2-in-1 Filter/O-Ring Insertion Tool Kit



2-in-1 Filter/O-Ring Insertion Tool Handle

Inserting an O-Ring (ASE 100/150/200/300/350 systems), using 2-in-1 Filter/O-Ring Insertion Tool Handle.



Place the O-ring over the tip of the tool.



Insert the tool into the center hole of the extraction cell cap.



Press the tool firmly inside the cap until the O-ring snaps into place.



Filter Insertion Attachments

Instrument	Compatible cell body size
ASE 100/150/300/350	10 mL, 34 mL, 66 mL, and 100 mL
ASE 200	5 mL, 11 mL, 22 mL, 33 mL

Questions about this or any other Restek product? Contact us or your local Restek representative (www.restek.com/contact-us).

Restek patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.) Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.

© 2019 Restek Corporation. All rights reserved. Printed in the U.S.A.

www.restek.com

#803-04-001 Rev. date: 01/19



RESTEK
Pure Chromatography